

Product Specification



The PSA 10.555 Calibration Kit is used for the calibration of the PSA 10.525 Sir Galahad. The System includes a calibration vessel, thermometer, and a set of gas tight syringes. Hg vapour is drawn from the vessel and injected into the PSA Sir Galahad system. The user then tells the software the temperature and volume of Hg vapour injected and the software will then calculates the Hg vapour injected based on the vapour pressure of the Hg.

The PSA 10.555 Calibration Kit is comprised of the following items:

Qty	Part No.	Description
1	G523V002	Calibration Vessel (assembled)
1	Z360S524	50µl Syringe
1	Z360S531	250µl Syringe
1	Z360S533	1ml Syringe
1	Z360S535	5ml Syringe
1	Z550T002	Thermometer: -0.5/+40°C x 0.1 Grads (UKAS certificate)
1 pk	Z36N052	Sideport Needles (3 per pack)
1 pk	Z012S001	Septa: Silicone/PTFE 0.075in. x 12mm (144 per pack)

www.psanalytical.com

P S Analytical

